



SUNPOWER

CryoTel™ Family  
Free-Piston Stirling Cryocoolers

OPERATING INSTRUCTIONS

Model : CryoTel MT  
Version 4  
Issued: March 25, 2010

---

## Intellectual Property Declaration



Manufactured by Sunpower, Inc., under the following patents: U.S. Pats 4,583,364; 4,602,174; 4,623,808; 4,649,283; 4,805,408; 4,864,232; 4,866,378; 4,912,409; 4,926,123; 5,003,777; 5,148,066; 5,342,176; 5,385,021; 5,450,521; 5,457,956; 5,461,859; 5,496,153; 5,502,968; 5,525,845; 5,537,820; 5,592,073; 5,642,088; 5,642,622; 5,715,693; 5,749,226; 5,775,273; 5,873,246; 5,941,079; 6,035,637; 6,038,874; 6,170,442; 6,199,381; 6,293,184; 6,446,336 B1; 6,536,326; 6,684,637; RE 38,337; 6,782,700; U.S. patents pending. Also: Australia Patents 676,805; 677,518; 680,770; 685,997; 701,785; 709,315; 753,580; 754032. Brazil PI 950368-5; PI 960 8885-0; PI 9710742-5; PI 9713840-1; PI 9713840-1. Canada patent 2,184,473. EPO 0655120 1B issued in France, Italy, UK, Netherlands, Sweden and Germany (as DE 69329862.6-08). EPO 0693160 issued in Italy, UK, Netherlands France, Germany (as DE 69403468 T2); EPO 0754364 1B issued in France, Italy, UK, Netherlands and Germany (as DE 69526217.3-08). EPO 0783618 1B issued in France, Italy, UK, Netherlands and Germany (as DE 69518926.3-08). EPO 0878014 1B issued in France, Italy, UK, Netherlands and Germany (as DE 69611387.2-08). EPO 0885413 1B issued in France, Italy, UK, Netherlands and Germany (as DE 69627894.4-08). Italy 1297082. UK Patent No. 0,218,682; UK 2,334,307; UK 2,330,651. DBP No. 0,218,682 (Germany).

India Patents 177477, 178274, 185034; 185035. Republic of Korea patents 0202290; 0292453; 0301548; 0309486; 0320093. Mexico patents 184451; 194065; 197407; 201368. New Zealand Patents No. 263331; 282959; 319499; 302849; 517329. Singapore Patent P-No. 48360; 51842; 87422; 87423; 87424. Taiwan patent NI-77875.

Japan Patents 特許第1986438号 特許第3100163号 特許第3224232号 特許第3413658号

Patents pending in United Kingdom, Germany, Italy, France, Netherlands, Sweden, Brazil, Japan, Republic of Korea, India, China, Singapore.

---

Sunpower, the Sunpower graphic identifier, and CryoTel are trademarks of Sunpower, Inc. and are registered in the U.S. and other countries. Other products, logos, and Company names mentioned herein may be trademarks of their respective owners.

---

---

## Table of Contents

I.	<b>Introduction</b> .....	5
II.	<b>Precautions (Dos and Don'ts)</b> .....	6
III.	<b>Unpacking</b> .....	7
IV.	<b>Mechanical Mounting</b> .....	8
V.	<b>Attachment of Heat Load to Cold Finger</b> .....	8
VI.	<b>Cold End Vacuum</b> .....	9
VII.	<b>Feedthroughs</b> .....	10
VIII.	<b>Temperature Sensor</b> .....	11
IX.	<b>Heat Rejection</b> .....	13
	A. Water Jacket.....	13
	B. Copper Fins.....	14
	C. Conducting Solid.....	14
X.	<b>CryoTel Operation</b> .....	15
	A. Preparation.....	15
	B. Beginning Operation.....	15
	C. Operation Shutdown.....	17
XI.	<b>LED and 5 Pin Connector Summary</b> .....	18
XII.	<b>CryoTel Serial Communications Command Reference</b> .....	20
	A. Serial Interface Specifications.....	20
	B. Controller Connector Wiring Information.....	20
	C. CryoTel MT Terminal Emulator Commands.....	20
	D. Remote On-Off control.....	21
	E. Allowable Number Formats.....	22

---

## FIGURES

Figure 1: CryoTel MT .....	5
Figure 2: CryoTel Cryocooler.....	7
Figure 3: Pressure Vessel Mounting Holes.....	8
Figure 4: Weld Flange holes .....	8
Figure 5: Cold Tip M3 Mounting Holes.....	9
Figure 6: Power Cable.....	10
Figure 7: Rear Pressure Vessel End Plate.....	11
Figure 8: Temperature Sensor and Wiring Harness.....	11
Figure 9: Test Cap Assembly .....	12
Figure 10: Controller Temperature Sensor Connector.....	13
Figure 11: Water Jacket Heat Rejection.....	14
Figure 12: Cooling Fin Heat Rejection.....	14
Figure 13: CryoTel MT Control Board .....	16
Figure 14: Controller LED's and 5 pin Connector.....	18
Figure 15: Optocoupler Circuit.....	18

---

**This page intentionally left blank.**

---

## I. Introduction

Sunpower, Inc. is pleased to present this manual with your newly-purchased CryoTel model cryocooler. As the world leader in free-piston Stirling cryocooling, Sunpower fully supports its products and is pleased to hear of any comments or concerns from new or existing clients.

### CryoTel™ MT Operating Instructions

The Sunpower cryocooler is a precision instrument; robust in many ways but fragile in others. This operating manual provides standard-use instructions and precautions. Please post where technicians will be working with the unit. Please do not assume that if a warning or instruction is not included here, a particular test or application of this unit is acceptable. It may not be. If you have any questions about this unit, or about any tests or applications you intend to perform, please contact Sunpower support service at phone (740)-594-2221; email [cryo-support@sunpower.com](mailto:cryo-support@sunpower.com)



**CryoTel MT dimensions:** length with balancer: 223 mm (8.9 in); length without balancer: 183 mm (7.2 in); diameter: 73 mm (2.9 in);  
mass: 1.95 kg

**Figure 1:** CryoTel MT

---

## II. Precautions (Dos and Don'ts)

Although the CryoTel MT is mechanically robust, it can be damaged if not handled properly when in operation or when removing it from its packaging. The handling and operational **Dos** and **Don'ts** are listed below. Refer to the diagram of the CryoTel MT in Figure 2.


**CRITICAL! – Do not operate the cryocooler without adequate cooling at the heat rejection site. Overheating the cryocooler will cause permanent damage.**

### **DON'T**

1. Pick up the cryocooler by the cold finger
2. Set the cryocooler on the cold tip.
3. Allow the cold finger to be dented. The slightest dent will render the unit inoperable.
4. Drill holes, or in any other way puncture, or attempt to modify, the pressure vessel.
5. Operate the unit without proper cooling. Heat must be removed from the copper heat rejection area of the cryocooler. If the cooler is provided with external cooling fins, air must be forced over the fins and the flow path should not be obstructed.
6. Puncture or otherwise damage the copper service tube.
7. Subject the electrical feedthroughs to mechanical stress; i.e. axial or radial movements, axial loads, blows, or the like.
8. Mount the cooler by suspending it from the balance absorber mounting bolt.
9. Apply clamping pressure to the pressure vessel (labeled in Figure 2).
10. Rigidly attach the absorber stud to “ground.” Instead, let the cooler “float” via a rubber bushing or other means of articulation.
11. Remove the protective cover on the cold weld on the end of the copper charge tube. Do not subject the cover to blows.
12. Control power to the cryocooler by making or breaking the power leads between the controller and the cryocooler.
13. Use an external, automatic, closed loop control system which attempts to control the cryocooler operation by varying the setpoint settings in the cryocooler controller.

### **DO**

1. Check with Sunpower before making any modifications to the cryocooler
2. Note that if the control board is installed in an enclosure, cooling must be provided.
3. If the cooler is operating in a laboratory test mode with heaters providing the thermal load to the cold tip, interlocks should be provided so that the heater cannot operate unless the cryocooler is running. This will prevent accidental overheating of the cold tip.

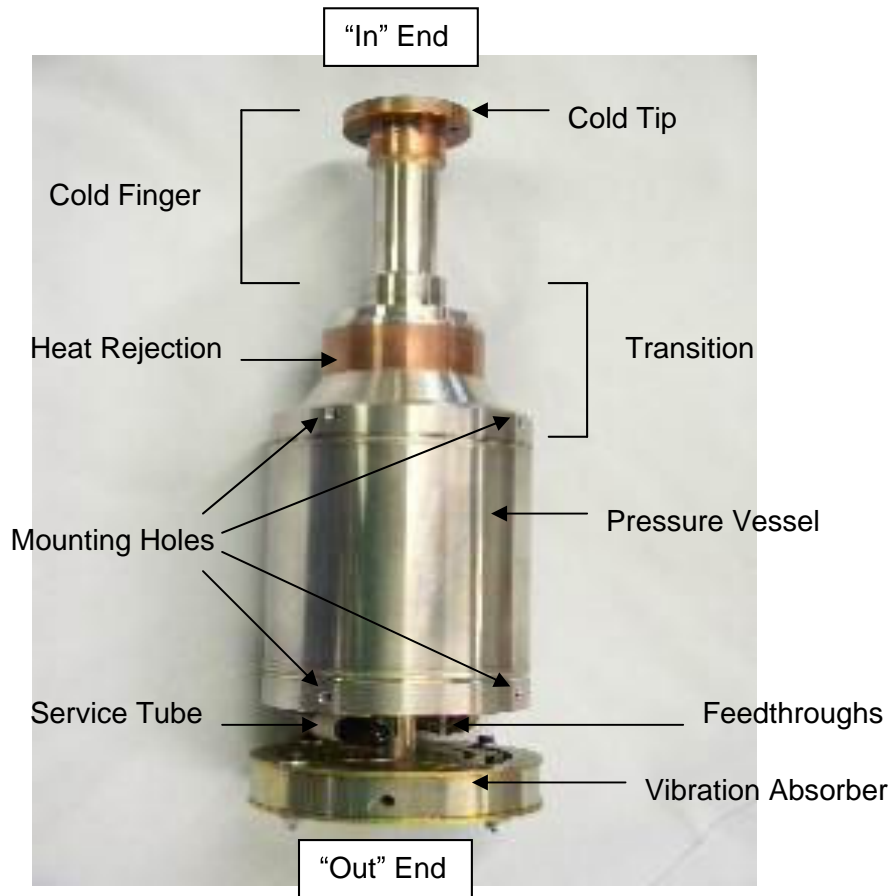

**Figure 2:** CryoTel™ Cryocooler

### III. Unpacking


When unpacking or handling, only hold the cryocooler by the pressure vessel or the transition (see Figure 2). Do not hold or support the cryocooler by holding the cold finger. The CryoTel™ MT may sit horizontally on a level bench top or table, but the cooler must be supported to prohibit rolling. The cold finger must be protected from any contact as well. The cold finger is delicate. Any deformation of the cold finger, such as small dents, bending, etc. will ruin the cryocooler.

---


## IV. Mechanical Mounting

The CryoTel™ MT has a thick metal end plate on the “out” end of the pressure vessel nearest the passive vibration absorber. M4 threaded holes on the outer diameter of the transition and end plate are designed for attachment of the cooler to an external mechanical structure for testing or integration into the application. These can be seen in Figure 3.

On the face of the external copper ring that rejects heat from the cooler is a stainless steel plate. The plate has four M3 threaded holes that can be used for attaching a removable NW50 or customer specified vacuum flange (see Figure 4). A vacuum flange can also be welded to this plate, but must be done with a low energy weld such as plasma or micro TIG. It is recommended that, if needed, the customer allow Sunpower to perform the welding operation. The mechanical structure including the cryostat attachment plate, heat rejection copper, and transition have been designed to allow mounting the entire assembly by the cryostat and cantilevering the cooler and vibration absorber from the cryostat in a given application.

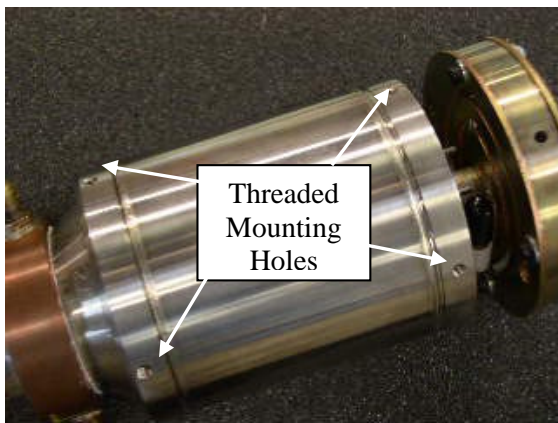

**Figure 3:** Pressure Vessel Mounting Holes

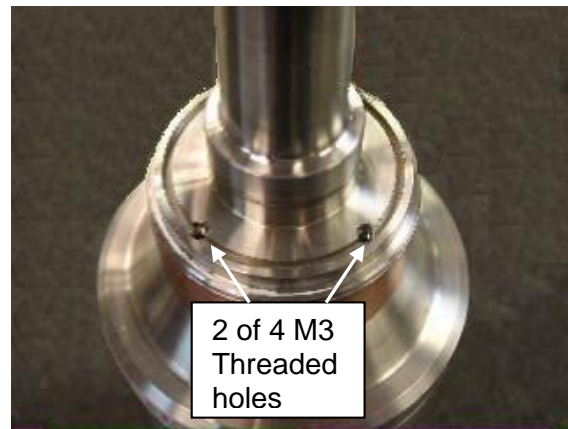

**Figure 4:** Weld Flange Holes

## V. Attachment of Heat Load to Cold Finger

The copper cold tip at the end of the cold finger has four M3 threaded holes that can be used for mounting equipment to the cold tip for applications or tests (see Figure 5). The application of force to the cold tip must be handled with care and in accordance with these not-to-exceed limits:

- Cantilevered force, cooler **not** in operation: do not exceed a cantilevered force on the cold finger (static or shock) > 100 N (10.20kg).
- Cantilevered force, cooler **in** operation: do not permit the heat load apparatus to deliver force to the cold finger (static or shock) > 30 N (3.05 kg)
- Vertical loading: do not exceed a pure tension or compression load > 300 N (30.50 kg)
- Torque, cold finger: do not exceed a load >16 N\*m (142 in\*lbs)



**Figure 5:** Cold Tip M3 Threaded Holes

## **VI. Cold End Vacuum**

The cold finger of the CryoTel™ MT operates best in an insulated vacuum. The vacuum eliminates possible loading of the cold finger from convection or condensation of elements in the atmosphere such as water vapor and nitrogen. The vacuum is created by a customer provided vacuum Dewar or cryostat. As noted above, a flange can connect the cryostat and the stainless steel vacuum flange with an o-ring seal or welded attachment (Fig 2).

When using a mechanical vacuum pump it is appropriate to seal the flange with an o-ring and clamp. The vacuum pump will remove any residual gasses released by the o-ring and continue to maintain a vacuum inside the cryostat. However, an o-ring may corrode and become brittle over an extended period of time.

For customer applications requiring long-term vacuum without using a vacuum pump, it is recommended that the cryostat be welded to the CryoTel vacuum interface.

---

## VII. Feedthroughs

The electrical pins on the metal plate at the end of the pressure vessel near the balance absorber are surrounded by glass, which acts as an insulator and prevents helium from leaking out of the pressure vessel. This arrangement is called a feedthrough. Because of the glass insulator, the feedthroughs should be handled cautiously. They are a permanent feature of the pressure vessel back end plate and should not be modified in any way.

The CryoTel MT is shipped with a power cable that attaches to the feedthroughs and the controller. This wiring harness consists of orange and white 16 gauge wires with a molded plastic connector for the feedthroughs at one end and a plastic connector housing that plugs directly into the controller output cable of the same wire colors. This harness is shown in Figure 6.



**Figure 6:** Power Cable


It is necessary to remove the balance absorber in order to install the power cable. The balance absorber is attached with the center M5 screw. Remove this screw and the balance absorber will be free from the cryocooler.

Install the power cable by aligning the cryocooler feed through pins with the holes in the cable connector and pressing down. Insert the retaining screw into the connector and tighten.

Install the balance absorber by re-installing the M5 screw.

### **Important assembly notes:**

- The heads of the four screws that fasten the balance absorber assembly together should be facing towards the cryocooler. The nuts for these screws will be on the “top” of the balance absorber.
- The screw heads of the four balance absorber assembly screws need to be positioned so that none of them are directly over either the power cord connector or the service tube elbow. This is to ensure that the screw head does not impact either of these features during large amplitude displacements of the balance absorber.

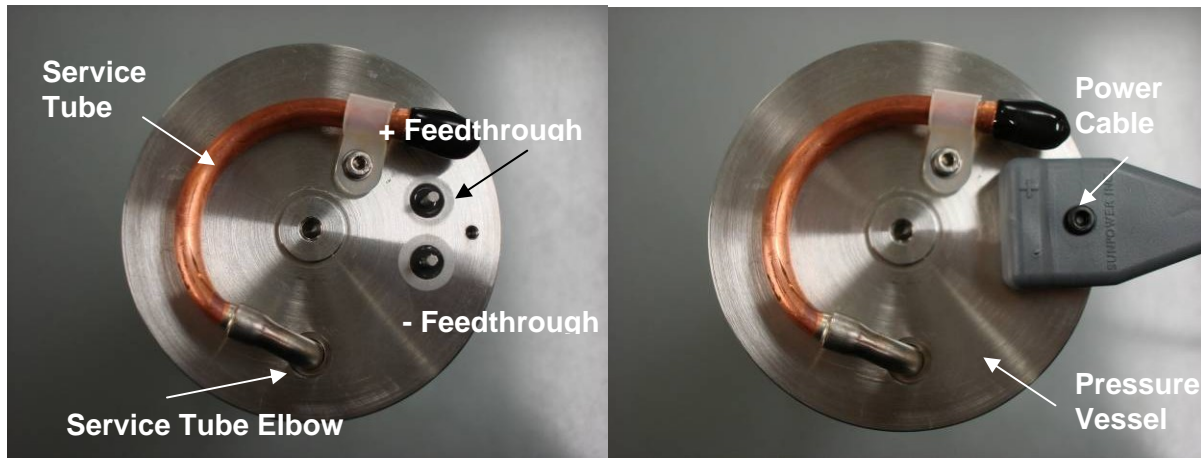


**Figure 7:** Rear Pressure Vessel End Plate and Power Cable Wire Harness

### VIII. Temperature Sensor

For proper operation of the CryoTel™ controller it is necessary to mount a temperature sensor to the object being cooled. The sensor allows the controller to measure and control the cooled object's temperature. The sensor feedback also controls the cooler's power ramp-up. If the sensor is not installed properly, the cooling capacity of the cryocooler will be severely limited and temperature control will not function.

The controller sensor may be either a PT-111 platinum RTD or equivalent. Sunpower packages the PT-111 in a copper disk so that the sensors are identical in size, shape and wiring. Figure 8 shows the sensor and the wiring harness connecting sensor to controller.

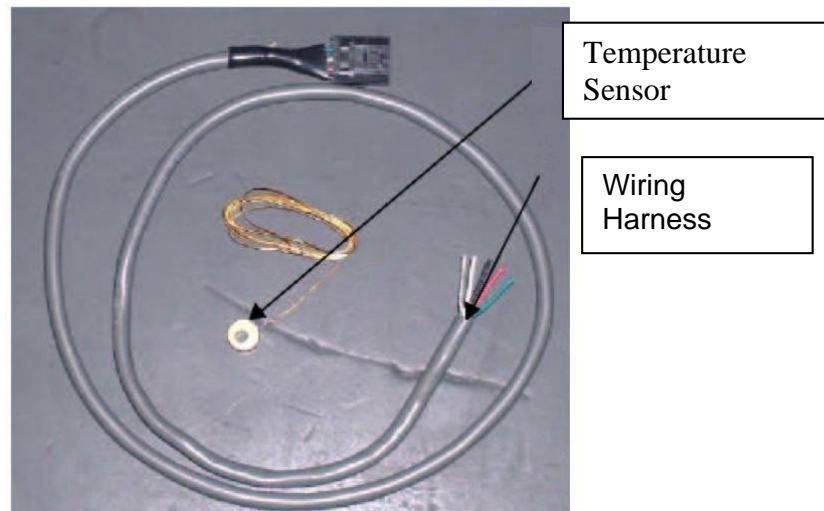


**Figure 8:** Temperature Sensor and Wiring Harness

The temperature sensor is encased in a small copper slug. Use an M3 screw in the 3mm drilled hole to attach the sensor to the test cap or to the object being cooled. Use a thin layer of

---

Indium, Apiezon grease, or similar thermal grease between the sensor and the object to insure proper thermal conduction.

Laboratory testing can be accomplished with a test cap arrangement similar to the one shown in Fig 9. The test cap consists of a copper disk, two resistors, two through holes with M3 clearance for connection to copper tip, and two M3 holes for attachment of sensors to the test cap. Calculate the heat load (power in Watts = current x voltage) by providing a known voltage and current across the resistors on the test cap. Mount the control sensor using one of the M3 mounting holes on the test cap. If using a second sensor, mount it to the test cap for independent temperature measurement. Apply an appropriate thermal grease to the bottom face of the test cap in order to maximize conduction. Mount the test cap to the cold tip using the M3 threaded holes on the cold tip.

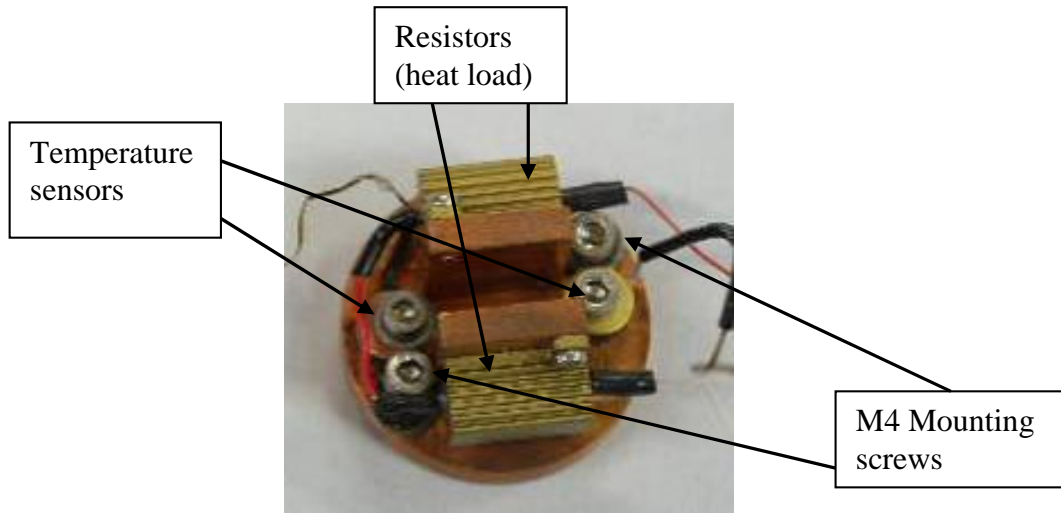


**Figure 9:** Test Cap Assembly

When the cold tip is contained in a vacuum vessel, use feedthroughs to connect the wiring harness to the sensor wires. Use at least four conductors connecting the feedthrough to the control temperature sensor. Use additional feedthroughs if employing a test cap or other sensors inside the cryostat. All wires inside the cryostat should be as small as possible in order to minimize parasitic loading on the cooler. The wires of the sensor and wiring harness are color coded to help insure proper connection. Table 1 shows the proper connection of the wires. Figure 10 and Table 2 describe proper connections of the wiring harness to the CryoTel Controller.

RTD Wire Color	Wiring Harness Color
Black	Black
Red	Red
Clear	White
Green	Green

**Table 1**

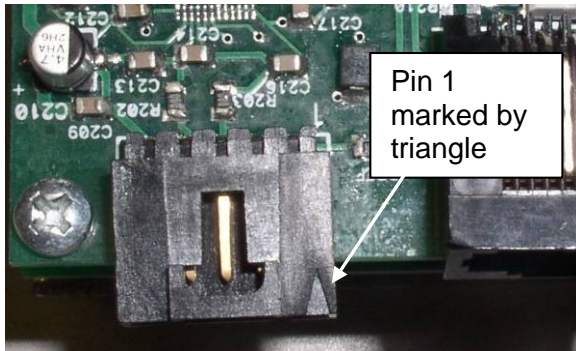

**Figure 10:** Controller Temperature Sensor Connector

Mating connector for temperature interface cable is DigiKey part # WM2903.

Terminals are DigiKey part # WM2554.

PIN	Function	RTD Wire Color	Harness Wire Color
1	I+	White (Clear)	White
2	I-	Red	Red
3	V+	Black	Black
4	V-	Green	Green
5	Shield		

**Table 2**

## IX. Heat Rejection

Proper heat rejection is crucial to the operation of a cryocooler. Some CryoTel heat rejection options are a water jacket, copper fins, or a conducting solid.

**CRITICAL!** – Do not operate the cryocooler without adequate cooling at the heat rejection site. Overheating the cryocooler will cause permanent damage.

### **A. Water Jacket**


Sunpower can provide a permanent or removable water cooling jacket with the CryoTel MT (Fig. 11 –both are identical in appearance). Use a suitable thermal grease to maximize conduction between the heat rejector and the water jacket if using a removable unit (the permanent unit is installed at the factory and does not require thermal grease). The thermal grease should be refreshed periodically as it will dry out over time. After connecting the water supply to the water jacket for heat rejection, ensure that air is removed from the water jacket. This can be done by flowing water through the cooling system for several minutes prior to the run. If it appears that air is still trapped in the water jacket, try carefully tilting the cooler in the direction away from the cooler discharge tube. This will raise the level of the discharge tube and allow any trapped air to escape. Trapped air will decrease heat rejection and cooler performance. Water should be flowing through the jacket at approximately 5mL per second (0.08 gallons per minute).



**Figure 11:** Water Jacket Heat Rejection

### **B. Copper Fins**

Copper fins are also available in a permanent or removable form. If copper fins are used for heat rejection (Fig. 12-*permanent fins are pictured*), use a fan with a rating of 100 cubic feet per minute ( $1.6\text{e-}3 \text{ m}^3/\text{min}$ ) for air through-flow. Sunpower recommends mounting the fan to a cylindrical shroud that can be placed around the fins in order to direct the air through the fins. Note that some heat rejection also occurs from the walls of the pressure vessel. Therefore avoid permitting any containment structure for the cryocooler to hinder the heat rejection from any part of the cooler. It is also recommended to electrically interlock the cooling fan and the controller in order to ensure that the cooler cannot operate without also turning on the fan. For additional cost, Sunpower is able to provide a fan, shroud, and controller that are suitable for laboratory testing.

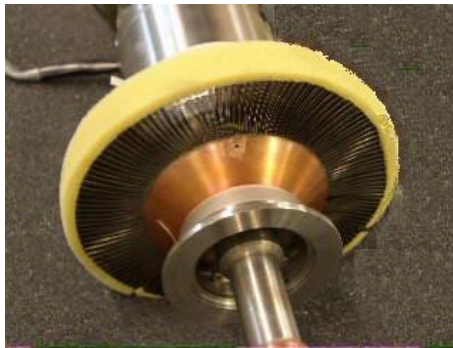

**Figure 12:** Cooling Fin Heat Rejection

### **C. Conducting Solid**

Use of any other means of heat rejection necessitates ensuring that the proper amount of heat is being transferred away from the rejector and that the reject temperature is suitable. It is recommended to involve Sunpower with the design of a conduction heat exchanger.

---

## X. CryoTel Operation

**CRITICAL!** – Do not operate the cryocooler without adequate cooling at the heat rejection site. Overheating the cryocooler will cause permanent damage.

### **A. Preparation**


1. Attach application apparatus or test cap to cold tip of cold finger.
  - a. Apply a thin layer of Apiezon thermal grease or Indium to the copper tip of the cold finger in order to maximize conductivity.
2. Attach the customer-provided cryostat by appropriate means if a vacuum is required for the application.
3. If using a vacuum, make wire connections from temperature sensor and any other measuring devices inside the cryostat through Dewar feedthroughs.
  - a. For maximum performance, wires should not touch the walls of the Dewar. Contact with walls will increase the heat loading (conduction) during use and will yield lower available cooling power.
4. Seal the cryostat vacuum except for the connection to the vacuum pump (if using vacuum).
5. Mechanically attach the cryocooler assembly to the customer provided mounting device.
6. Connect the wiring from the cryostat feedthroughs to the appropriate test and measurement devices.
  - a. The cold tip temperature sensor must be connected to the controller's mating connector as described previously using the provided wiring harness and connector (See Figure 13 for controller connections).
7. Attach the vacuum pump to cryostat, start the vacuum and allow it to reach a vacuum of  $10^{-4}$ Torr or better.

### **B. Beginning Operation**

1. Prepare the CryoTel MT for testing as described above.
2. Connect the power cables for the CryoTel MT to the controller's mating connector (controller connections in Fig. 13).
  - a. Verify the correct polarity is observed with respect to the CryoTel MT – orange power cable wires are positive, white wires are negative.
  - b. Attach the included wire harness to the cooler as shown in Figure 6 then attach the plastic connector at the other end to the connector with orange and white wires on the controller (Fig 13).
    - i. Note: The connector attached to the cooler wires is an AMP Mate-N-Lok® plug housing, 1-480698-0; with AMP Mate-N-Lok® contacts, 3505471-1 (available from Mouser Electronics, part numbers 571-14806980 and 571-3505471 respectively).
    - ii. Note: The connector attached to the controller wires is an AMP Mate-N-Lok® cap housing, 1-480699-0 with AMP Mate-N-Lok® contacts, 3505501 (available from Mouser Electronics, part numbers 571-14806990 and 571-3505501 respectively).
3. Connect the controller to a 24V<sub>DC</sub> power outlet.
  - a. Note: The input power connector attached to the control board is an AMP Mate-N-Lok® cap housing, 1-480699-0 with AMP Mate-N-Lok® contacts, 350547-1 (

available from Mouser Electronics, part numbers 571-14806990 and 571-3505501 respectively).

- b. Note: Use a suitable connector for the power supply such as the same AMP Mate-N-Lok® plug housing used for the cryocooler power cables, AMP part number 1-480698-0 with AMP universal connector socket, 350550-1.
  - c. The controller has a startup sequence that positions the piston in the axial center. This sequence takes approximately 7-10 seconds once the controller is turned on.
  - d. The controller then applies a 60 Hz AC voltage to the motor, creating the oscillating motion of the piston that begins the cooling process.
  - e. The controller increases the power input to the cryocooler based on cold end temperature and predetermined time constants. The maximum draw by the controller is approximately 100 W<sub>E</sub>.
4. Allow cold tip temperature to stabilize at desired cold end temperature
    - a. Proper operation of the CryoTel MT requires a minimum loading of 2 Watts. If minimum load requirements are not met, the stabilized temperature will be below 77 K.

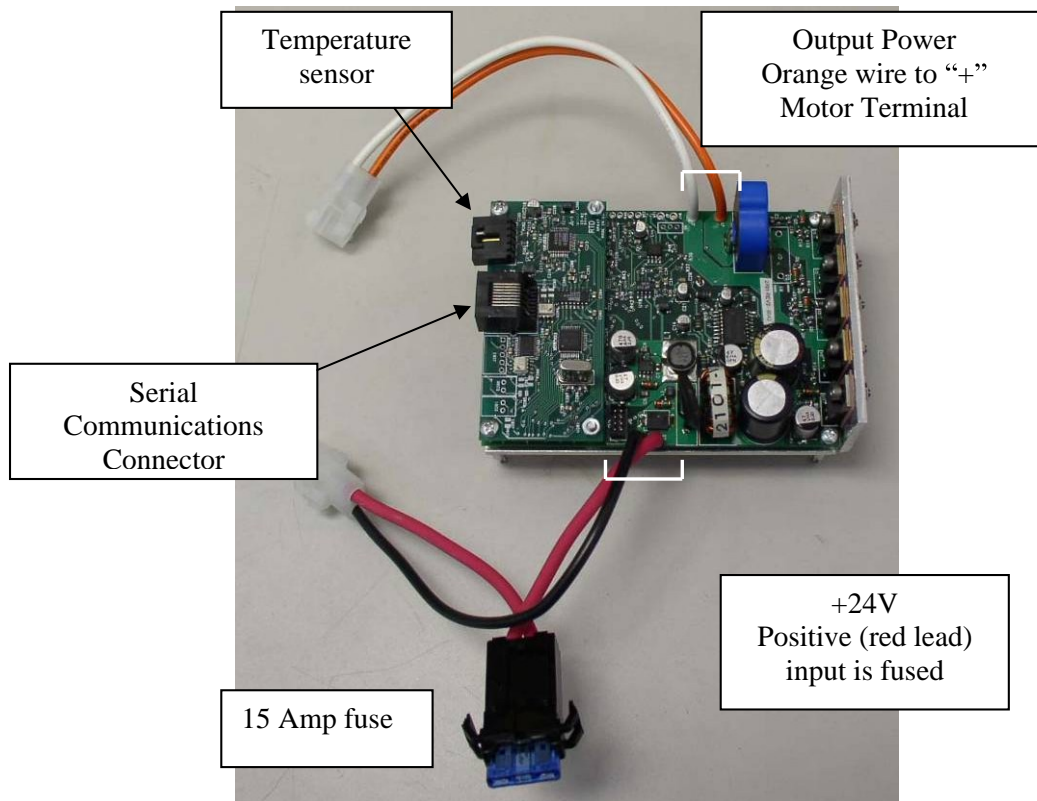

**Figure 13: CryoTel MT Control Board**

---

### **C. Operation Shutdown**


1. Turn off the controller.
  - a. The quick rattle heard when power is cut off from the CryoTel is due to the stored energy in the vibration absorber being dissipated.
  - b. In rare cases where the operating temperature was below 60 K, it has been observed that the cooler can energize itself during shutdown and become an engine. To prevent this, allow the cold tip to reach a temperature above 60 K before shutting down.
2. Allow the cold tip and the rest of the cryocooler temperature to rise to room temperature before opening the vacuum in order to prevent the vapor outside the cryostat from condensing and freezing on the cold tip.
  - a. Applying a heat load to the cold tip during the warming process reduces the time needed to reach room temperature.
  - b. Sunpower recommends using a relay in conjunction with the temperature sensor to ensure that the power supply to the heat load shuts off when the cold tip temperature reaches 300 K.

---


## XI. LED and 5 Pin Connector Summary

The two LED's, shown in figure 13, indicate whether the cooler has reached its set point temperature or if the cooler is still in cool down mode. The red LED will be on as long as the cooler is in cool down mode. When the cooler reaches its set point temperature within 0.5K the green LED will turn on.

The 5-pin connector is connected to an optocoupler circuit as shown in Figure 15. When the green LED is on, current is able to flow from pin 1 to pin 5 through the Darlington Phototransistor. When the green LED is off, the current flow through the phototransistor is blocked. This would be useful if the user wanted switch an externally powered indicator light or control system input on when the cooler reaches the set point temperature. Pins 2, 3, and 4 on the 5 pin connector are not used.

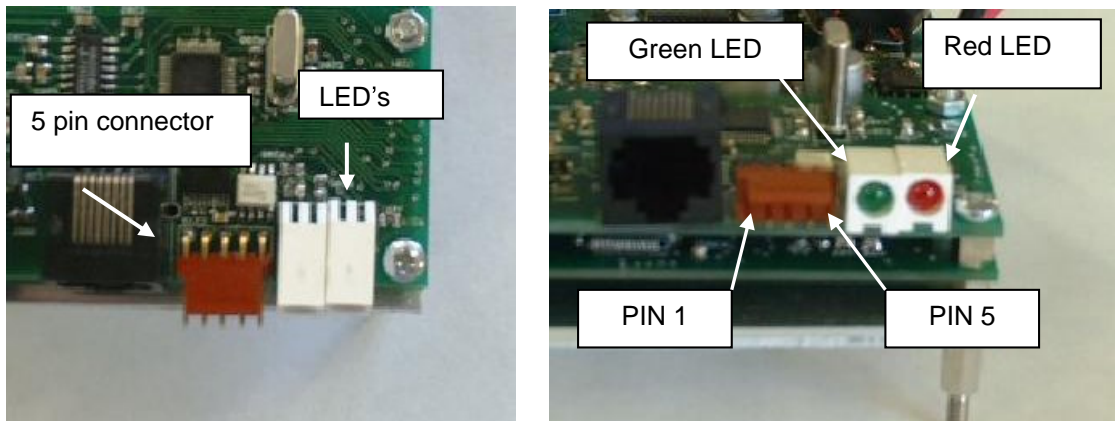

**Figure 14:** Controller LED's and 5 pin Connector

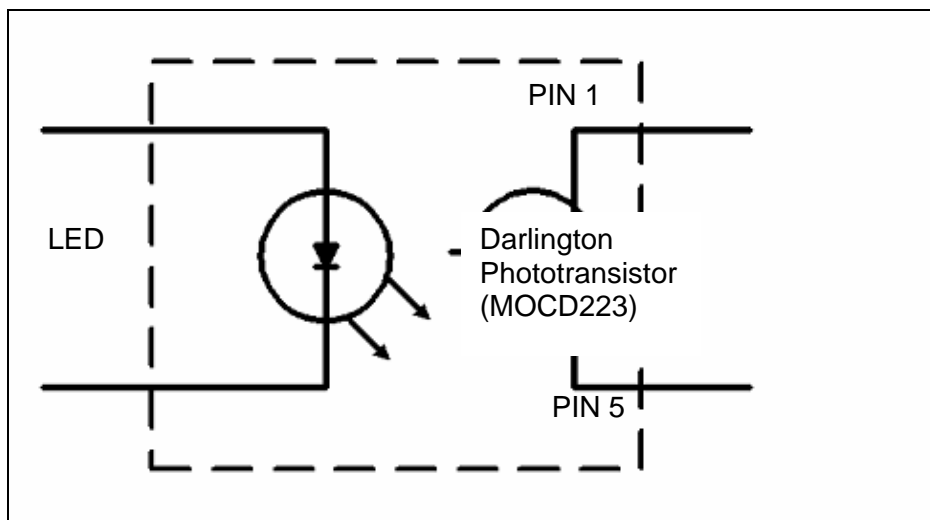

**Figure 15:** Optocoupler Circuit.  
Phototransistor is on when green LED is on.

---

**This page intentionally left blank.**

---

## XII. CryoTel Serial Communications Command Reference

Control the CryoTel™ MT with any terminal emulator program. A common choice is Hyperterminal, as it is included with Windows. The connecting cable and computer adapter are included.

**Caution:** Use of an external program, such as LabVIEW™, to establish closed loop control of the cryocooler via serial communication, may cause controller malfunction. This malfunction condition requires the controller to be returned to Sunpower, Inc. for reprogramming. Manual setpoint adjustments may be made as required without problem.


### A. Serial Interface Specifications

RS-232  
Baud Rate: 4800  
Flow control: none  
Data bits: 8  
Stop bits: 1

### B. Controller Connector Wiring Information

Pin Numbering of RJ-45 connector and Pin-out of RS-232 Connector on CryoTel Controller

Pin	Function
1	RX
2	GND
3	Do not connect
4	GND
5	Do not connect
6	GND
7	TX
8	GND

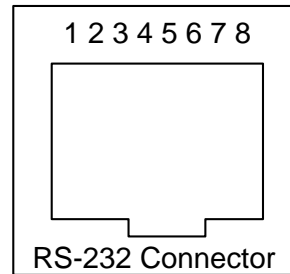

### C. CryoTel MT Terminal Emulator Commands

The following pages show an explanation of each of the terminal emulator commands and display examples.

1. Display temperature sensor value in Kelvin
  - a. Command: **TC<CR>** (<CR> means “press Enter key”)

```
TC
330.00
```

6. Show the current controller mode
  - a. Command: **SET PID<CR>**
  - b. The value returned will be a two or a zero. If it is a zero, the controller will maintain a constant input power as commanded by **SET PWOUT**. If it is a two, the controller will try to maintain a constant temperature as commanded by **SET TTARGET**.

```
SET PID
002.00
```

7. Set controller Mode
  - a. Command: **SET PID=V<CR>**
  - b. V may be a 2 or a 0 as shown before. 2 is for closed-loop temperature control and a 0 is for manual, input power control.
  - c. When this command is entered, the return display is equivalent to using SET PID without a specific value.

---

```
SET PID=0
000.00
SET PID=2
002.00
```

4. Display target temperature in Kelvin
  - d. Command: **SET TTARGET<CR>**
  - e. This command shows the target temperature of the controller when in mode 2

```
SET TTARGET
077.00
```

5. Set target temperature in Kelvin when in mode 2
  - a. Command: **SET TTARGET=V<CR>**
  - b. V corresponds to the target temperature. This value is stored in flash so it is retained in the controller at power up.

```
SET TTARGET=86.42
086.42
```

6. Display command stroke when in mode 0 control.
  - a. Command: **SET PWOUT<CR>**

```
SET PWOUT
150.00
```

7. Set command power when in mode 0 control
  - a. Command: **SET PWOUT=V<CR>**
  - b. V corresponds to the target power in watts. While any number from 0.0 to 999.99 can be input, the controller will only command a power that will not damage the cryocooler.
  - c. The minimum power is 30W @ 77 K.
  - d. The maximum power is a function of cold head temperature and increases as the cold temperature decreases.

```
SET PWOUT=80
80.00
```

8. Display current command power and power limits
  - a. Command: **E<CR>**
  - b. The top value is the maximum allowable power for the current temperature. The middle value is the minimum allowable power. The bottom is the current command power. All values displayed with this command are in watts.

```
E
090.00
030.00
080.00
```

9. Remote ON-OFF feature:

There is a control feature that is available that allows the power output to the cryocooler to be disabled while keeping the controller powered up. The controllers are shipped with this function disabled. The majority of our customers does not need it and would not want to make the wiring changes to connect the pins. Pins 3 and 5 on the RJ45 serial communications port are used as the signal input. Shorting these pins together enables the controller output; open circuit disables the controller output.

---

You can enable this feature by sending these commands to the controller:

Unlock <enter>  
SSTOP=1 <enter>  
Lock <enter>

<<Caution>> It is imperative that the “Lock” command be entered at the end of the sequence to prevent the software being compromised during controller operation.

Note: after you enable this function, pins 3 and 5 must to be connected in order to drive the cryocooler.

**D. Allowable Number Formats**

When using the terminal emulator to command the controller, all input numbers should be in one of the following formats. X is a digit from 0 to 9.

X	XX	XXX
X.X	XX.X	XXX.X
X.XX	XX.XX	XXX.XX

---

**End of CryoTel® MT Operating Instructions Version 4**